



Starting-up your Zymocube

- 1 Turn on the Zymocube
- 2 Start the Zymosoft
- 3 Scan QR code
- 4 Type your experiment name
- 5 Open the tray
- 6 Place your zymplate on the tray
- 7 *Optional: Activate the advanced mode*
- 8 Click on Next

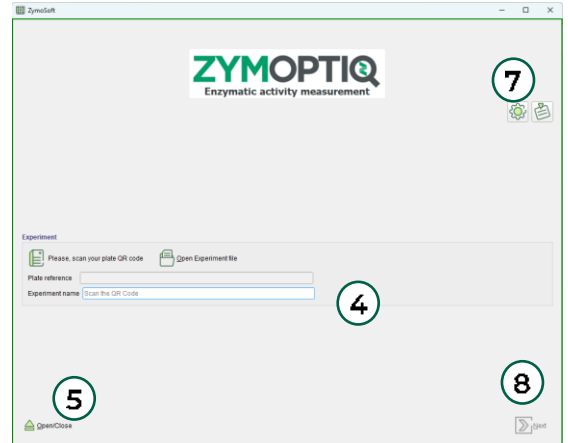

Acquiring your data

- 1 Add your experimental data (create it, edit a template or upload a file)
- 2 Define well types, areas, add information
- 3 Click ok
- 4 Launch readout

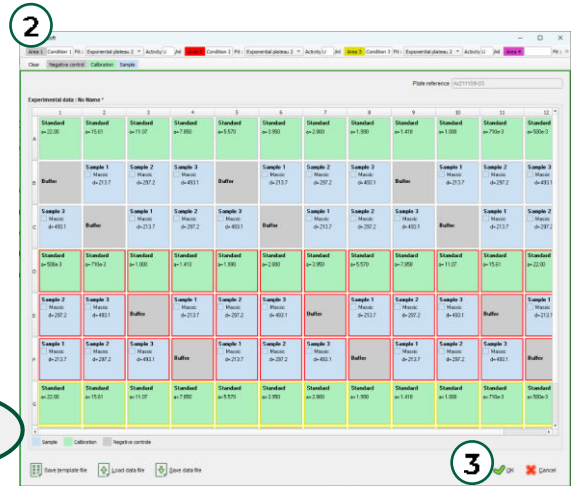

Analyzing your results

Check your results:

- 1 Negative control
 - 2 Calibration curve and data
 - 3 Results per sample
 - 4 Data details per well
- (you can navigate between the different views)
- 5 Repeat for all areas
 - 6 Open tray to retrieve your plate
 - 7 End your Experiment



Data is automatically saved